

PW21 The Analysis of Alprazolam in Synthetic Urine via Dual Column Switching

Charles Yang, Jonathan Beck and Rohan Thakur
Thermo Electron Corporation, San Jose, CA.

Alprazolam, Column, Mass Spec, SPE, Online sample prep, benzodiazepine

Analysis of urine for Alprazolam without sample preparation using 2 column switching system automated by a software program.

The quantification of Alprazolam as a drug of abuse in urine has traditionally been performed by GC/MS, followed SPE extraction. This requires substantial time in sample preparation and run time. A novel method utilizing 2 columns, a switching valve, 2 LC pumps, and a triple quadrupole mass spectrometer is described here. This method requires no sample clean up prior to injection onto the LC column.

Injection of synthetic human urine containing Alprazolam and Alprazolam-d5 were injected onto 2 Thermo Hypersil GOLD reverse phase column. Two pumps were used; one to elute the compound, the second to re-equilibrate and wash the off line column. Upon elution of the analytes, the valve is actuated and the next injection is performed. A triple quadrupole mass spectrometer, Thermo Scientific TSQ Quantum Discovery MAX, is used to analyze compounds.

A standard curve was prepared in synthetic human urine (UriSub, CST Technologies, Inc.) covering ranges 7.8 to 500ng/mL. Three quality controls were prepared at 11.7, 187.5 and 750ng/mL and a second set of standards was prepared at 1:10 dilution of the calibrators and QCs using di-water. The calibration curves are linear over both of the ranges described above. The columns showed no deterioration in quality or performance. Even at 1:10 dilution, the calibration curve is linear and well above the detection limit of the instrument.